



Now offering online ordering (USA only) of many popular products through PayPal secure shopping.

Reference Photos of Phototubes:

[View Common Phototube and C-mount Adapters.](#)

| | | |
|--|--|--|
|  <p>28mm ID</p> <p>American Optical Series 10 28mm ID Photoport without #1044 Phototube. MDSLR-AO fits this port directly.</p> |  <p>23mm ID 25mm OD</p> <p>AO 10 showing #1044 Phototube. Older #1044's are usually silver. MM25TC clamp recommended for MM Series adapters. SC67LJ or SC50LJ C-mount adapters recommended for dedicated microscope cameras.</p> |  <p>23mm ID (25mm OD)</p> <p>Meiji ML Series has standard 23mm ID / 25mm OD phototube. MM25TC clamp recommended for MM Series adapters. SC67LJ or SC50LJ C-mount adapters recommended for dedicated microscope cameras.</p> |
|  <p>39mm</p> <p>Leica HC dovetail Photoport - usually marked "HC" and uses a 3mm allen screw. Uses DDxxHCC C-mounts. MDSLR-HC fits this port directly.</p> |  <p>37mm ID</p> <p>Leica DMR with 37mm ID Photoport</p> <p>Looks very similar to the 38mm type, so measure carefully! Uses D10DMC or DD63MZC or DD50MZC C-mounts. MDSLR-DMR can be special ordered to replace this phototube for DSLR's.</p> |  <p>36mm</p> <p>Leica DME 36mm ID Photoport. MDSLR-LE fits this port directly.</p> |
|  <p>38mm ID</p> <p>Older Leitz ISO 38mm ID Photoport - used for most Leitz trinocular scopes: Ortholux, Orthoplan, Dialux, Laborlux, etc. Uses DxxNLC 1x C-mount for 38mm ISO port. Uses DD20NLT ISO T-mount. MDSLR-LM fits this port directly.</p> |  <p>Leica M and S series stereomicroscopes with externally threaded port with 28mm ID. Uses Leica C-mounts or Leica 10445930 conversion to 37mm ID phototube plus DDxxMZC C-mounts.</p> <p>MDSLR-LM fits this port directly.</p> |  <p>37mm</p> <p>Leica MZ 37mm ID Phototube HU/HV</p> <p>Off-white phototubes are 37mm. Uses D10DMC or DD63MZC or DD50MZC C-mounts.</p> |
|  <p>38mm ID</p> <p>Wild M3 with ISO 38mm ID Phototube</p> <p>Black Leica/Wild Phototubes are 38mm. Uses DxxNLC C-mounts. Uses DD20NLT ISO T-mount.</p> |  <p>30mm ID</p> <p>Older Wild M5 with 30mm Phototube - older ones are painted yellow</p> <p>Direct fit for MM Series; dedicated microscope cameras require SCxx30 C-mounts.</p> | |

| | | |
|---|---|--|
|  <p>Nikon Eclipse 55i with Ergo Head and built-in 0.7x C-mount Phototube. Works with 1/2" and 2/3" sensor cameras directly, but no way to attach DSLR or MM series adapters. Also may restrict some C-mount cameras due to size constraints.</p> |  <p>Nikon Eclipse LV100 with 40mm ID Photoport. Requires Nikon MBB63430 to convert to standard 38mm ISO port for use with DxxNLC C-mounts. MDSLR-NK fits this port directly. This is the new Nikon standard port for materials and stereo microscopes.</p> |  <p>Nikon E200 with non-ISO 38mm port. Requires Nikon MBB73520 Y-T TV Tube for use with DxxNLC C-mounts. MDSLR-NE fits this port directly.</p> |
|  <p>Nikon Eclipse with ISO 38mm ID Photoport - Nikon "chimney" tube has been removed. Uses DxxNLC C-mounts. Uses DD20NLT ISO T-mount.</p> |  <p>Nikon Eclipse with "Chimney" phototube with recessed 30mm ID port. MM-SLR, MM99 and MMC00L and variants fit this chimney tube directly; MM3 does not. Uses SCxxNF C-mounts.</p> |  <p>Nikon Labophot-2 ISO 38mm Photoport with "chimney" tube removed. Uses DxxNLC C-mounts. Uses DD20NLT ISO T-mount.</p> |
|  <p>Nikon Optiphot / Labophot "chimney" type phototube with recessed 23mm ID port. MM-SLR, MM99 and MMC00L and variants fit this chimney tube directly; MM3 does not. Uses SCxxNF C-mounts.</p> |  <p>Olympus BH2 with reversed dovetail photoport with recessed 23mm ID port. For MM Series, use MMBH2 clamp. Uses DxxBHC C-mounts.</p> |  <p>Olympus BX with 51mm ID dovetail Photoport. Uses DxxBXC C-mounts. MDSLR-BX fits this port directly. Also for Olympus SZX.</p> |

| | | |
|--|---|--|
|  <p>27mm ID Martin Microscope Company</p> <p>B&L Stereozoom 7 with 27mm ID photoport. Older SZ7's accept a threaded 25mm phototube.</p> |  <p>28mm ID Martin Microscope Company</p> <p>B&L / Leica Stereozoom 6 with 28mm ID photoport. Contact us for help with this one.</p> |  <p>48mm dovetail Martin Microscope Company</p> <p>Jenaval 48mm dovetail Photoport. Jenalum and Jenavert also use this port. Uses DxxZJC C-mounts. Jenapol's also have this port, but the phototube must be much shorter for parfocality, so it requires M10ZIP C-mount - 1x only.</p> |
|  <p>30mm ID Martin Microscope Company</p> <p>Zeiss Stemi 2000C with new 30mm ID photoport (Zeiss # 45 29 02). Uses DxxZNC C-mount or MDSLR-ZS DSLR Adapter.</p> |  <p>Zeiss Axio Imager side photoport uses newer Zeiss DxxZNC C-mounts or MDSLR-ZS DSLR Adapter.</p> |  <p>Zeiss Axioskop with 44mm dovetail photoport. Uses DxxZAC C-mounts.</p> |
|  <p>Zeiss Axioplan with 44mm dovetail photoport. Uses DxxZAC C-mounts.</p> |  <p>47mm ID Martin Microscope Company</p> <p>Zeiss Universal 47mm dovetail Photoport. Uses DxxZUC C-mounts.</p> |  <p>23mm ID 25mm OD Martin Microscope Company</p> <p>Zeiss Standard with 23mm ID / 25mm OD phototube. For MM Series, MM25TC clamp recommended. Uses SCxxLJ C-mounts.</p> |

[Home](#) - [Microscopes](#) - [Brands](#) - [Digital Imaging](#) - [Company History](#) - [Contact Information](#)

Copyright 2001- 2013 Martin Microscope Company All rights reserved.
Prices, specifications, and availability subject to change.